

Milli-Q Gradient and Milli Q Synthesys (1996-2002)

The cabinet of the system needs to be removed. If the system is wall mounted, unhook the system from its wall mounting.

1. Put the System in STANDBY mode by pressing on OPERATE/STANDBY for 2 seconds.
2. Switch of power of the system by putting the main swifch (to position 0)
- 3 Unplug and disconnect the main elecfrical power cord of the System
4. Open fhe Poinf-Qf-Use 3 way valve to depressurize the System.
5. Milli-Q Synthesis oniy: Unscrew and remove the cleaning porf plug (B).
6. Unscrew the two screws on the lower clips (C).
7. Slide the two clips backwards (D) to unlock the cabinet
8. Lift up the cabinet slighly, disconnect the ground wire and remove the cabinet
9. Remove the UV lamp power supply cord (E).
10. Using a screw driver, unscrew and remove the UV lamp housing head (F)
11. Use protective Gloves to replace the UV lamp.
12. Follow the above instrucfions in reverse order to reassemble the unit
13. In PRE-OPERATE mode, open the Poinf-Of-Use valve for 5 mn to rinse the unit

The uv lamp contain mercury this should therefore be disposed of by the owner

B) UV lamp lifetirne reset

In STANDBY or PRE-OPERATE mode

Press "Measure + Cleaning + Menu" buttons together to enter the Ser vice MENU
Press 4 times on MENU button

The UV lamp timer is reset for 365 days

MENU
start position.

Press on MENU buton for 2 seconds to return to OPERATE

The replacement procedure is now completed. The Milli-Q is ready for use.